


Application Note

G β γ -GIRK1 Interaction Assay



Introduction

AlphaScreen™ Gβγ-GIRK1 Interaction is a fully homogeneous assay in 384-well format, designed to measure interaction between G protein βγ subunit and the carboxyl terminal domain of the GIRK1 potassium channel (named cGIRK1). Gβγ is biotinylated and binds to streptavidin-donor beads, while cGIRK1 binds to anti GST-acceptor beads via its GST tag. Specific interaction between the two proteins leads to signal increase. The binding is fully reversible and completely abolished by competition with G protein α subunit. This assay can be easily applied to any protein-protein interaction.

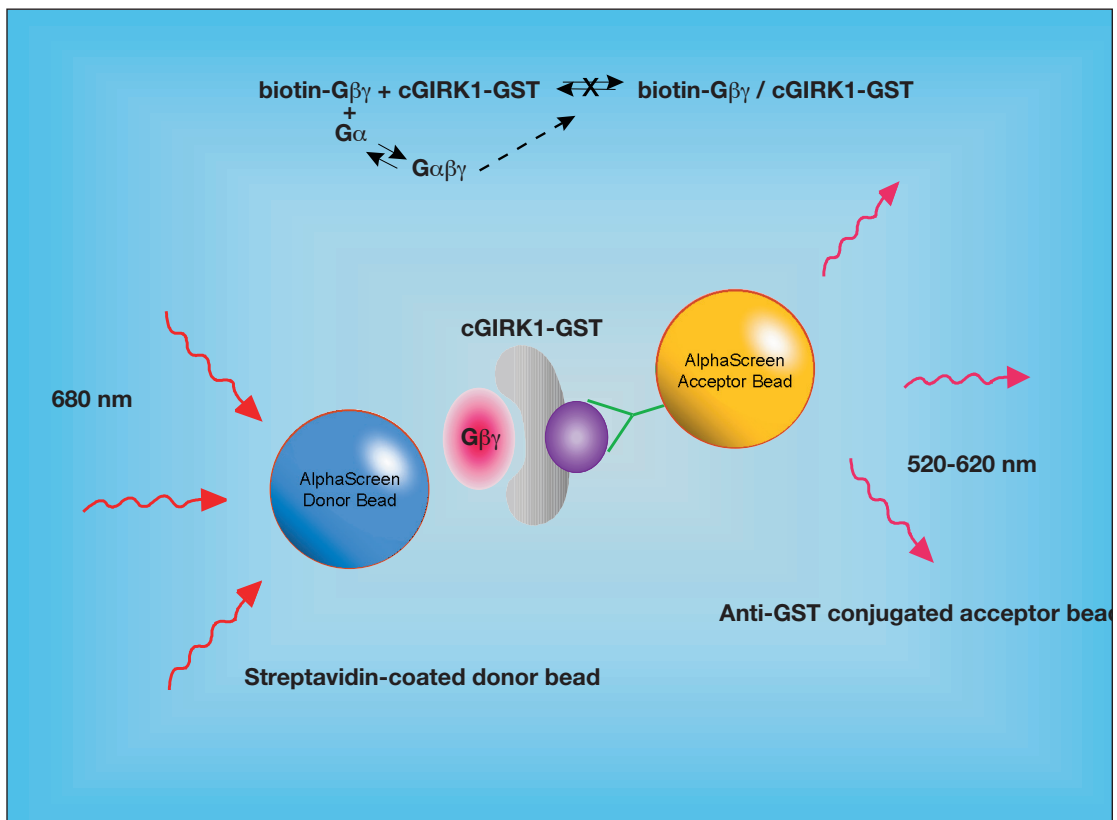

Example #1:


Competition of biotin-G β 1 γ 2 binding to GST-cGIRK1 by G β 1 γ 2

Assay developed in PerkinElmer 384-well OptiPlate™

Reagents

1. Biotinylated G β 1 γ 2 (12.3 μ M in 20 mM Hepes, 100 mM NaCl, 0.1% CHAPS): dilute to 125 nM with assay buffer
2. GST-cGIRK1 (14.7 μ M in 25 mM Hepes pH 8.0, 50 mM NaCl, 5% glycerol): dilute to 50 nM with assay buffer
3. G β 1 γ 2 (22.1 μ M in 20 mM Hepes, 100 mM NaCl, 0.7% CHAPS): dilute to 30 μ M - 5 nM with assay buffer
4. Streptavidin-donor beads (5 mg/ml in 25 mM Hepes, pH 7.4): dilute to 50 μ g/ml with assay buffer
5. Biotin-G β 1 γ 2 / streptavidin-donor bead complex: mix equal volumes of biotinylated G β 1 γ 2 and streptavidin-donor beads and incubate for 30 minutes at RT
6. Anti GST-acceptor beads (5 mg/ml in 25 mM Hepes, pH 7.4): dilute to 125 μ g/ml with assay buffer

Assay buffer: 100 mM Hepes, pH 7.4, 100 mM NaCl, 0.1% CHAPS, 1 mM EDTA, 0.1% BSA

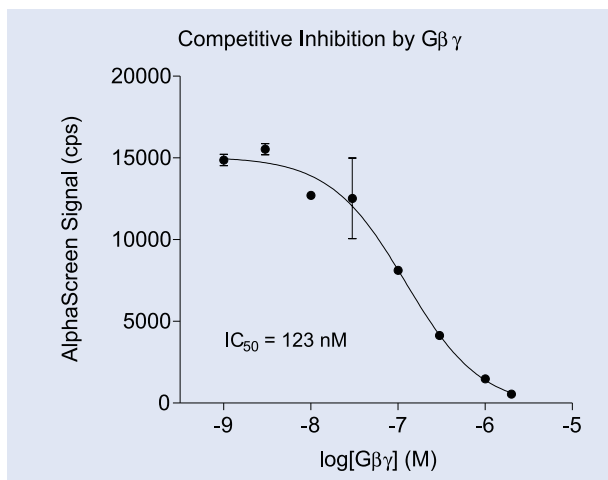

Protocol:

1. Add 5 μ l GST-cGIRK1.
2. Add 5 μ l G β 1 γ 2.
3. Add 5 μ l anti GST-acceptor beads.
4. Incubate 30 minutes at RT.
5. Add 10 μ l biotin-G β 1 γ 2 / streptavidin-donor bead complex.
6. Incubate 60 minutes at RT.
7. Read plate.

Example #2:


Inhibition of biotin-G β 1 γ 2 binding to GST-cGIRK1 by G α _i

Assay developed in PerkinElmer 384-well OptiPlate™

Reagents

1. Biotinylated G β 1 γ 2 (12.3 μ M in 20 mM Hepes, 100 mM NaCl, 0.1% CHAPS): dilute to 125 nM with assay buffer
2. GST-cGIRK1 (14.7 μ M in 25 mM Hepes pH 8.0, 50 mM NaCl, 5% glycerol): dilute to 50 nM with assay buffer
3. G α _i1 (11.5 μ M in 20 mM Hepes, 100 mM NaCl, 0.7% CHAPS): dilute to 3 μ M with assay buffer, incubate for 2h at 30°C in 25 mM Tris-HCl, pH 7.5, 0.5 mM MgCl₂, 100 μ M GDP, then further dilute to 2.5 μ M - 5 nM with assay buffer
4. Streptavidin-donor beads (5 mg/ml in 25 mM Hepes, pH 7.4): dilute to 50 μ g/ml with assay buffer
5. Anti GST-acceptor beads (5 mg/ml in 25 mM Hepes, pH 7.4): dilute to 125 μ g/ml with assay buffer
6. GST-cGIRK1 / anti GST-acceptor beads complex: mix equal volumes of GST-cGIRK1 and anti-GST-acceptor beads and incubate for 30 minutes at RT

Assay buffer: 100 mM Hepes, pH 7.4, 100 mM NaCl, 0.1% CHAPS, 1 mM EDTA, 0.1% BSA

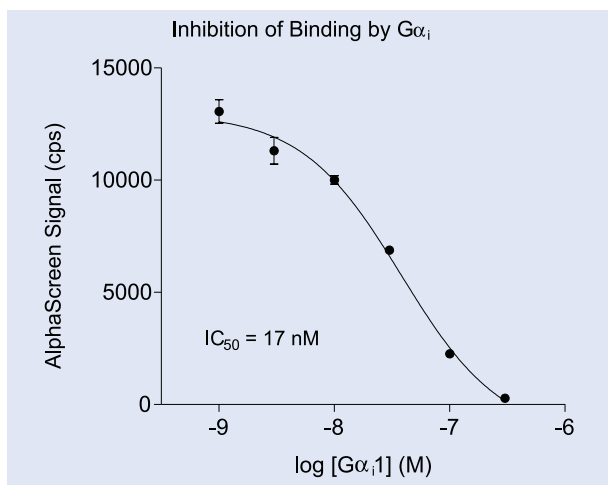

Protocol:

1. Add 5 μ l biotin-G β 1 γ 2.
2. Add 5 μ l G α _i1.
3. Add 5 μ l streptavidin-donor beads.
4. Incubate 30 minutes at RT.
5. Add 10 μ l GST-cGIRK1 / anti-GST-acceptor bead complex.
6. Incubate 60 minutes at RT.
7. Read plate.

Example #3:


Reversibility of biotin-G $\beta\gamma$ / GST-cGIRK1 interaction

Assay developed in PerkinElmer 384-well OptiPlate™

Reagents

1. Biotinylated G $\beta\gamma$ 2 (12.3 μ M in 20 mM Hepes, 100 mM NaCl, 0.1% CHAPS): dilute to 125 nM with assay buffer
2. GST-cGIRK1 (14.7 μ M in 25 mM Hepes pH 8.0, 50 mM NaCl, 5% glycerol): dilute to 50 nM with assay buffer
3. G $\beta\gamma$ 2 (22.1 μ M in 20 mM Hepes, 100 mM NaCl, 0.7% CHAPS): dilute to 5 μ M with assay buffer
4. Streptavidin-donor beads (5 mg/ml in 25 mM Hepes, pH 7.4): dilute to 50 μ g/ml with assay buffer
5. Anti GST-acceptor beads (5 mg/ml in 25 mM Hepes, pH 7.4): dilute to 125 μ g/ml with assay buffer

Assay buffer: 100 mM Hepes, pH 7.4, 100 mM NaCl, 0.1% CHAPS, 1 mM EDTA, 0.1% BSA

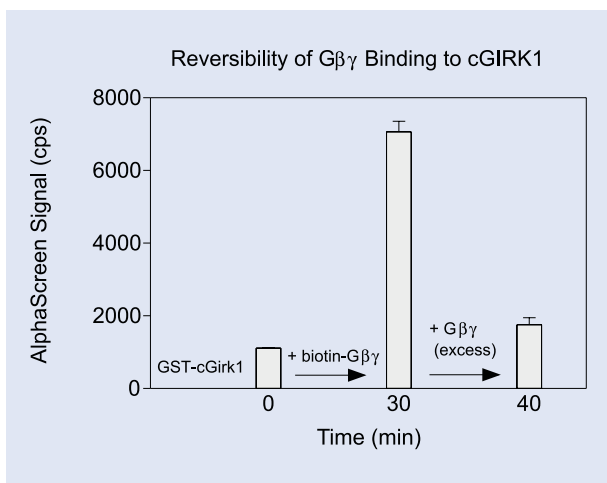

Protocol for binding association:

1. Add 5 μ l GST-cGIRK1.
2. Add 5 μ l anti GST-acceptor beads.
3. Incubate 30 minutes at RT.
4. Add 5 μ l biotin-G $\beta\gamma$.
5. Incubate 30 minutes at RT.
6. Add 10 μ l streptavidin-donor beads.
7. Incubate 10 minutes at RT.
8. Read plate.

Protocol for binding dissociation:

1. Add 5 μ l GST-cGIRK1.
2. Add 5 μ l biotin-G $\beta\gamma$.
3. Add 5 μ l anti-GST-acceptor beads.
4. Incubate 30 minutes at RT.
5. Add 5 μ l G $\beta\gamma$.
6. Incubate 10 minutes at RT.
7. Add 5 μ l streptavidin-donor beads.
8. Incubate 10 minutes at RT.
9. Read plate.



PerkinElmer Life and Analytical Sciences, 710 Bridgeport Avenue, Shelton, CT 06484 USA (800) 762-4000 or (+1) 203-925-4602

PerkinElmer Life and Analytical Sciences, Imperiastraat 8, BE-1930 Zaventem Belgium

Technical Support: in Europe: techsupport.europe@perkinelmer.com in US and Rest of World: techsupport@perkinelmer.com

Belgium: Tel: 0800 94 540 • **France:** Tel: 0800 90 77 62 • **Netherlands:** Tel: 0800 02 23 042 • **Germany:** Tel: 0800 1 81 00 32 • **United Kingdom:** Tel: 0800 89 60 46
Switzerland: Tel: 0800 55 50 27 • **Italy:** Tel: 800 79 03 10 • **Sweden:** Tel: 020 79 07 35 • **Norway:** Tel: 800 11 947 • **Denmark:** Tel: 80 88 3477 • **Spain:** Tel: 900 973 255

© 2003 PerkinElmer, Inc. All rights reserved. PerkinElmer is a registered trademark and AlphaScreen and OptiPlate are trademarks of PerkinElmer, Inc. PerkinElmer reserves the right to change this brochure at any time without notice and disclaims liability for editorial, pictorial, or typographical errors. Printed in USA.

006815 [M3337 6/00]

www.perkinelmer.com